

Stereo Microscope System



For Life Science Use

Comfortable, High-Quality Life Science Imaging





User Comfort Is Critical for Precision Performance

The SZX7 stereo microscope is easy to use and delivers outstanding optical performance so that users are comfortable performing imaging tasks from advanced research to routine inspections. The microscope's Galilean optical system, previously restricted to more specialized microscopes, offers high zoom ratio, as well as high image clarity, true-to-life color, and fidelity reproduction of the specimen in crisp, well-defined detail.

The SZX7 microscope can be customized using a range of accessories to accommodate a variety of specimen types and sizes.



Quality Optics for Consistently Superior Image Reproduction

View your specimens clearly and accurately without straining your eyes. With its smooth finish, apochromatic zoom optics, eyepieces, and easy adjustability, the SZX7 minimizes strain and fatigue while fulfilling the key mission of Olympus microscope designers—to provide the optimal image for any specimen. The performance of the high-level Galilean optical system is complemented by much less distortion than before with a high numerical aperture (NA).

The SZX7 microscope body is manufactured using lead-free optics, demonstrating Olympus' commitment to protect the environment.

7:1 Wide Zoom Ratio

With a magnification range of 8X–56X (using a 1X objective/10X eyepieces), the SZX7 microscope offers a zoom ratio of 7:1. This high zoom ratio enables a specimen to be observed at appropriate magnification.

Excellent Resolving Power

High-quality objectives deliver accurate, high-resolution images that show specimens in minute detail.

A Range of Objectives to Suit Every Specimen and Every Application

• Superior Image quality with high resolution and excellent flatness: The DFPLAPO1X-4 objective provides excellent optical performance with plan apochromat correction and an NA of 0.10.

• Long working distance (W.D.):

Objectives range from the SZX-ACH1X (90 mm W.D.) to the DFPL0.5X-4 (171 mm W.D.). As a result, difficult to access surfaces can be easily observed.

Accurate Color Reproduction

Careful selection of lens surface coatings and apochromatic zoom optics make it possible to observe and document specimens with accurate color reproducibility.



Observation tubes



Galilean optics feature two (right/left) independent and parallel zoom optical paths. This system enables high optical performance as well as system modularity.

Convenient Front-Access Operation

Improved access to the most frequently used knobs and controls maximizes operator comfort and reduces neck and back strain.

Quickly Recall Magnification Settings via Click-Stop Mechanism

Many inspection and documentation tasks require the use of a known zoom magnification setting to get consistent and comparable results. The integrated click-stop mechanism provides quick and easy access to this important function.



Objectives



Tilting trinocular tube



OLYMPUS

XL

Fluorescence Unit with Coaxial Illumination Enables Clear, Bright Observation Even with Weakly Emitting Specimens

This reflected light fluorescence unit is used to observe fluorescence in living cells under a stereo microscope. The high-performance fluorescence filter sets have sharp cutoffs and high transmission to capture even faint emissions from fluorescent proteins.

Reflected Light Fluorescence Unit SZX-RFA

Three fluorescence filter blocks can be mounted in a 4-position slider. An open position is provided for easy access to transmitted light observation. The light source is a 100 W mercury lamp, for bright fluorescence observation with high contrast. A total of six filter sets are available, depending on the purpose.

High-Performance Filters for GFP/YFP

Two different types of high-performance filter sets are available for GFP/YFP. Optimized for the characteristics of GFP/YFP wavelengths, they have high transmission rates of 90% to 95% and sharp cutoffs for efficient detection of even weak fluorescence.



Filter cubes















Choose the Illumination Source That Suits Your Sample

LED Illuminator Stand / SZ2-ILST

The LED stand features a thin design to keep sample positions low and to optimize usability. Simultaneous transmitted and reflected light are available on this stand. LED light offers both a long lifetime and consistent color temperature at any intensity.



Transmitted Illumination Attachment / SZ2-ILA

Used with the SZ2-ST this cost-effective illumination stand provides bright, uniform illumination from low to high magnifications. An adjustable mirror provides direct and oblique illumination for low contrast specimens. An available LED light source (SZ2-CLS or BX3M-LEDT) provides the necessary power for a variety of illumination needs.



Choice of Your Suitable and Observation Method / SZX2-ILLTQ/SZX2-ILLTS

With a slim 41.5 mm (1.6 in.) design that is approximately half the thickness of previous halogen lamp transmitted light illumination bases, the LED transmitted light illumination bases have a lower height to enable a low eyepoint and easy access to basemounted samples during observation and operation. The LED illumination base SZX2-ILLTQ with quad position turret enables the user to choose cartridges and to switch from brightfield (standard/ high/low), oblique (standard/high/low), darkfield, polarized illumination, and shutter with a simple turn. A one position LED illumination base is also an option (SZX2-ILLTS). This makes the SZX2 series a flexible all-in-one microscope for various samples and observation tasks. Another advantage of LED illumination is a cooler base surface, which is suitable for long duration manipulation of live specimens. Power consumption is lower than a conventional 30 W halogen light source. A lifetime of over 60,000 hours significantly reduces operation costs.





Product	Observation Methods and Contrasts		
 SZX2-CBFL 	Brightfield, low-contrast		
2 SZX2-CBF	Brightfield, standard		
③ SZX2-CBFH	Brightfield, high-contrast		
(4) SZX2-COBL	Oblique, low-contrast		
5 SZX2-COB	Oblique, standard		
6 SZX2-COBH	Oblique, high-contrast		
⑦ SZX2-CSH	Shade plate		
⑧ SZX2-CDF	Darkfield		
③ SZX2-CPO	Polarization plate		



A variety of fiber guide illumination systems are available.



Flexible light guide / SZ2-CLGSF



Dual interlock light guide / SZ2-CLGDI



Six-point ring light guide / SZ2-CLGR



Coaxial reflected light illuminator / SZX2-ILLC10

Flexibility from Digital Imaging to Observing Large Specimens



DP74 Digital Camera

The DP74 color fluorescence camera captures realistic, highquality images and has features that enable users to make their observations easily. With a wide field of view, operators can capture images of more of their sample, quickly.

Light Beam Splitter / SZX2-LBS

Two digital cameras can be attached simultaneously. The light path can be changed between three different settings: 100% observation, 100% digital camera, and 50% observation and 50% to both left and right cameras.

Photo Adapter / SZX-PHA

Various adapters are available for different kinds of CCD cameras and can be used with the beam splitter.



Ergonomic Tilting Trinocular Tube / SZX2-LTTR Extendable Eyepoint Adjuster / SZX2-EEPA

The tilting trinocular tube and eyepoint adjuster enable comfortable microscope work, even over long periods. Users can adjust the tilting tube angle from 5 to 45 degrees and move the height of the eyepoint within a 120 mm range to maximize comfort. Moreover, attaching a digital camera enables users to obtain high-resolution images at 1920 \times 1440 pixels, which exceeds standard high-definition resolution.



Side-by-Side Discussion Tube / SZX-SDO2

Ample distance (650 mm) is provided between the primary and secondary observers, making observations easy without disturbing microscope operation. The color of the built-in pointer can be changed to contrast with the specimen.



Various Universal Stands

A variety of universal stands are available for the observation of large specimens. No matter the size of your sample, Olympus has the right choice of stands to suit any requirements.



SZX7



SZX7+SZ2-ST



SZX7+SZ2-ILST





69

SZX7 (SZX2-TR30 configuration) +SZ2-ILST



SZX7 specifications

ltem		Specifications						
Zoom microscope body SZX-ZB7		Zoom drive: Horizontal knob system Click stop for each zoom magnification: ON-OFF switching possible Zoom ratio values: 7:1 (0.8X to 5.6X) Zoom magnification indication: 0.8, 1, 1.25, 1.6, 2, 2.5, 3.2, 4, 5, 5.6 Objective mounting: screw mounting into thread Lead-free						
		Aperture iris diaphragm control: T	he AS unit (SZX-AS) is mountable	9				
Observati	on tube	SZX-BI45	SZX2-TR30	SZX2-TR30PT	SZX2-LTTR*1			
SZX-BI45 SZX2-TR3 SZX2-TR3 SZX2-LTT	30 30PT	View inclination angle 45°	Filting binocular tube Trinocular tube Ergonomic Long Tilting T View inclination angle 30° View inclination angle 30° View tilting angle 5° to 45 Light path selection: 2 steps Light path selection: 2 steps Light path selection: 2 steps Binocular 100%, (Binocular 100%, (Binocular 50%/Photo 50%)					
		All observation tubes: Lead-free		111010 10070				
	Interpupillary distance adjustable range	52 to 76 mm Eyepiece clamping knob provideo	1	57 to 80 mm Eyepiece clamping knob provider				
Extendable	e Eyepoint adjuster	SZX2-EEPA: Height adjustment ra	ange: 30–150mm, (with a scale a	ttached)				
Stand SZ2-ST SZ2-ILST		SZ2-ST	SZ2-ILST					
		Standard stand	LED reflected/transmitted illumination stand					
	Frame installation	Mounting diameter 76 mm						
	Focusing adjustment	Knob rotation tension adjustment Focusing stroke 120 mm						
	Stage plate	SZ2-SPBW (Black & white) SP-C (Glass clear transparent)	100 mm diameter dedicated glass plate is included					
	Light source	Compact light guide illuminator (SZ2-CLS) mountable (option) Transmitted light illumination attachment (SZ2-ILA) mountable (option)	Transmitted illumination: LED Reflected illumination: LED Average LED life span: 6000 hrs. Input rating: 100–120 V/200–240 V~0.15/0.1 A, 50/60 Hz					
Objectives	s	Model		NA	Working distance			
		DFPL0.5X-4*2 DFPL0.75X-4 DFPLAPO1X-4 SZX-ACH1X DFPLAPO1.25X-2 SZX-ACH1.25X DFPL1.5X-4 DFPL2X-4 All objectives: Lead-free		0.05 0.075 0.10 0.10 0.125 0.125 0.125 0.15 0.20	171 mm 116 mm 81 mm 90 mm 60 mm 68 mm 45.5 mm 33.5 mm			
Eyepieces	6	WHSZ series All eyepieces: Lead-free						
Weight	Configuration 1	4,360 g (9.6 lb)	5,200	g (11.5 lb)	5,300 g (11.7 lb)			
				6,000 g (13.2 lb) 6,100 g (13.4 lb)				

*1 SZX2-LTTR: intermediate magnification is 1.25X
 *2 The SZ2-ET auxiliary sleeve is required when the SZ2-ST/SZ2-ILST is used
 Configuration 1: SZX-ZB7 + DFPLAPO1X-4 + individual observation tube + WHSZ10X-H (2) + SZ2-ST
 Configuration 2: SZX-ZB7 + DFPLAPO1X-4 + individual observation tube + WHSZ10X-H (2) + SZ2-ILST

■ SZX7 optical performance*³

Eyepiece	epiece WHSZ10X-H WHSZ10X		WHSZ15X-H		WHSZ20X-H WHSZ20X		WHSZ30X-H	
FN	22		16		12.5		7	
Objective	Total magnification	Field of view (mm)	Total magnification	Field of view (mm)	Total magnification	Field of view (mm)	Total magnification	Field of view (mm)
0.5X	4X-28X	55–7.8	6X-42X	40.0–5.7	8X–56X	31.3–4.5	12X-84X	17.5–2.5
0.75X	6X-42X	36.7–5.2	9X-63X	26.7–3.8	12X–84X	20.8–3.0	18X–126X	11.7–1.7
1X	8X-56X	27.5–3.9	12X-84X	20.0–2.9	16X–112X	15.6–2.2	24X–168X	8.8–1.3
1.25X	10X-70X	22-3.1	15X–105X	16.0–2.3	20X-140X	12.5–1.8	30X-210X	7.0–1.0
1.5X	12X-84X	18.3–2.6	18X–126X	13.3–1.9	24X–168X	10.4–1.5	36X-252X	5.8–0.83
2X	16X-112X	13.8–1.9	24X–168X	10.0–1.4	32X-224X	7.8–1.1	48X-336X	4.4-0.63

*3 SZX2-LTTR: Intermediate magnification is 1.25X SZX2-ILLC10: Intermediate magnification is 1.5X

■ WHSZ eyepiece

	FN	Diopter adjustment	Reticle	Focal magnification
WHSZ10X	22	_	NA	—
WHSZ20X	12.5	_	NA	—
WHSZ10X-H	22	-8-+5	Yes*4	—
WHSZ15X-H	16	-8-+5	Yes*4	—
WHSZ20X-H	12.5	-8-+5	Yes*4	1.3X
WHSZ30X-H	7	-8-+5	Yes*4	2X

*4Applicable reticle size: 24 mm diameter, t1.5

SZX7 System Diagram



*1 Focusing unit (SZX2-FOF, SZX-FOFH or SZX-FO) and SZX-STF are required when mounting fluorescent unit.
 *2 SZX2-FOFH and SZX2-STL2 are required when using SZX-SDO2.
 *3 Please contact your nearest Olympus dealer for applicable cameras.
 *4 SZH-P400 and SZH-P600 can be attached to the transmitted light Illuminators.
 *5 Equipped to SZX2-ILLC10.



*6 Please contact your nearest Olympus dealer for applicable combination. *7 Not available in some areas. *8 SZX-ACH1.25X and DFFLAPO 1.00-** Common Section 270 II ST and S72-ST cannot be combined with SZX-RFA. *10 SZX2-STADM cannot be combined with SP-FL, SZ2-SPBW, SP-C and SZX-PO. *8 SZX-ACH1.25X and DFPLAPO1.5X-4 cannot be combined with SZX-STF.

Images are courtesy of National Institute for Basic Biology, Spectrography and Bioimaging Facility, Joe Sakamoto Ph.D., Yasuhiro Kamei Ph.D. (cover page, top right)

- OLYMPUS CORPORATION is IS014001 certified.
- OLYMPUS CORPORATION is ISO9001 certified.
- OLYMPUS CORPORATION is IS013485 certified.
- Illumination devices for microscope have suggested lifetimes.
 Periodic inspections are required. Please visit our website for details.

All company and product names are registered trademarks and/or trademarks of their respective owners.
 Images on the PC monitors are simulated.
 Specifications and appearances are subject to change without any notice or obligation on the part of the manufacturer.





OLYMPUS CORPORATION Shinjuku Monolith, 2-3-1 Nishi-Shinjuku, Shinjuku-ku, Tokyo 163-0914, Japan